Probe Software

Software for MicroAnalysis

Probe for EPMA

Probe Image

PictureSnapApp

Probe Software, Inc. 885 Crest Drive Eugene, OR 97405 USA (541) 343-3400 sales@probesoftware.com www.probesoftware.com



Correction for Secondary Boundary Fluorescence from Nearby Phases

Probe Software announces a new feature for the correction of secondary fluorescence boundary effects from nearby phases containing the element of interest that is detected by WDS or EDS spectrometers.

This boundary fluorescence effect is generally of concern for trace element analysis, but at distances less than tens of microns from phase boundaries, it can be a percent level artifact for some element couples, such as Cu-Co, Ni-Fe, Cr-W, etc.

Correction of these boundary fluorescence effects is possible through recent efforts by Francesco Salvat and Xavier Llovet at the University of Barcelona, Spain.

Starting with the development of sophisticated monte-carlo models for electron-photon fluorescence effects,



followed by new work implementing these insights into analytical expressions, means that these once time consuming and tedious operations can now be performed in a fraction of the time previously required for high precision quantitative calculations.

Probe Software has implemented these newly developed boundary fluorescence calculations in our CalcZAF acquisition and analysis software for the automatic correction of boundary effects in real-time based on the actual spatial distances and analyzed compositions.

	Most meny		Help Tal	ke-off In Kilos	rolts 15	0K
Select Material A	Select Material B	Select Material B Std		140	10	
(beam incident material)	(boundary material)	[primary standard]	DENIDARE Down			
2Magnani m malal	DDE Dauce metal	EVE David metal	DENERMA Prompt	Pennuor Panal B	Sinary Calculations	Close
13 Aluminum metal	505 Cabon (graphite)	505 Cabon (graphite)	* renermariant	Titemps Titemps		
14 Silicon metal	512 Magneoium metal	512 Magnesium metal	- Calculate Secondary Fb	assessment Profiles For the	Specified Element an	nd X-ray
21 Scandium metal	513 Aluminum metal	513 Aluminum metal		Design of the province of	0.00.0	
22 Titanium metal	514 Secon metal 521 Scandium metal	514 Secon metal 521 Scandium metal	Parameter	Files are created in PENEPHA,	Rootventiuor	
24 Chronium metal	522 Titanium metal	522 Titanium metal	Matenal A (beam incid	ent) Duper		Brow
25 Manganese metal	523 Vanadium metal	523 Vanadium metal	Material B (boundary)	Coper		Brow
25 iron metal	524 Diromium metal	524 Chromium metal				
22 Nickel metal	525 Manganese metal	525 Manganese metal	Material 8 Std (primary	coper		810
29 Copper metal	527 Cobalt metal	 527 Cobst metal 	Element Co y X-l	Ray ka y Microns 50	Points 50	Cos
Density Ose Strength Ose Econom	Density One Strength One 5	and Dentily Out Sharesth Out Fo	1	·	1	Clip
Denity Otc. Stenghotc. Energy	Denay Otc. Stenghotc. E	nergy Density Usc. Stenginusc. En	Bun Fanal (generate	k-ratio Exce	el 🔽 Use Grid Lines	_
96 U 10	lea lo lo	18'a lo lo	file for couple bou	ndary) 🛛 🗆 Use Log Scale	* Zoon Ful	
Oscillator strength (Fcb) a	and Oscillator energy (W/cb) of the pla	ismon should be zero for insulators				-
Create PENEPMA Material A	Create PENEPMA Materia	B Create PENEPMA Mat. B St	ad b			
From List	From List	From List	Co ka in	Cu adjacent to Co	(15 keV Cost	d)
				eu adjacem to co	(10 he 1, 00 h	-
Create PENEPMA Material A From	Cieste PENEPMA Material B Fr	om Create PENEPMA Material B Std Fr	hom _			
Formula	Formula	Formula			Cu (8.96) <	> Co (8.90
with dV anul: 5.8	Deputs of Sc equals 3	. Deschud Sc equile 3	K-Ratio %		-	
ensity of Crequels 7.19	Density of Tileguals 4.5	Density of Ti equals 4.5	Calc Wr %	(leah)	•	- 4.5
ensity of Mn equals 7.43	Density of V equals 5.8	Density of V equals 5.8	Baundan M	Di dicen Met Di	-	
ensity of Fe equals 7.86	Density of Crequals 7.19	Density of Cr equals 7.19	Boundary W	n. % (from Mat B)	-	-40
entity of Lo equals 0.9	Density of Mn equals 7.43 Density of Fe equals 7.85	Density of Mn equals 7.43 Density of Fe equals 7.85	 CalcZAP Wt 	.% (Armstrong/Love Sco	ott (derauit))	
maty of full equals 0.5	Density of Co equals 8.9	 Density of Co equals 8.9 	-			
					-	-3,5
					•	
	the second state of the second state of the	TATE OF A CONTRACT OF A DECIDENT OF A DECIDENT	CC		-	
inary Intensity Calculations (cro	pate .PAB files for one or all .M	IAT files) ("10 hours each at 3600 se				-3.0
inary Intensity Calculations (cre Create Material F	sate .PAB files for one or all .k les in PENEPMA_Path	(AT files) ("10 hours each at 3500 st Run Penlluor and Etail fi	w 1		•	- 3.0
inary Intensity Calculations (cm Create Material F 8. A Current	eate .PAB files for one or all .b lies in PENEPMA_Path	(AT files) ("10 hours each at 3600 sr Run Penlluor and Fital fi Browse ALL three materials	loe			-3.0
inary Intensity Colculations (cre Create Material P 8. A Current	eate .PAB files for one or all .b lies in PENEPMA_Path	(AT files) ("10 hours each at 3500 sr Bowse Bowse Current Bowse Current	lor			-3.0
Intensity Calculations (cm Create Material P R. A [Curred R. B [Correct	eate .PAB files for one or all .b lies in PENEPMA_Path	(AT files) (~10 hours each at 3600 sr Browse Browse				-3.0
Intensity Colculations (con- Create Material F 8. A Current 8. B Corrent 8. B Std Corrent	eate .PAB files for one or all .b iles in PENEPMA_Path	AT files) ("10 hours each at 3500 sr Browse				- 3.0 - 2.5 - 2.0
Internity Colculations (con Create Material P R. A [Current R. B [Corrent R. B Std [Corrent]	eate .PAB files for one or all .b	(AT files) (~10 hours each at 3500 sr Bowre ALL bree moterials Bowre Bowre Ban Perfluor/Filal for Materia Only Control Filal for Materia				- 3.0 - 2.5 - 2.0
Inary Intensity Colculations (cm Create Material F 8. A [Current 8. B [Cornel 8. B 514] [Cornel Time (n seconds per simul	ale .PAR files for one or all .N	AT files) (*10 hours each at 3500 st Bowre	100			- 3.0 - 2.5 - 2.0
Internity Calculations (cre Create Material F 2. A [Current 8. B [Correat [Current] 1. B Std [Correat Time (in seconds per simul Number of Simulations (in s	ate .x10	(AT files) (~10 hours each at 3500 sr Bowre Bowre Bowre Bowre Pan Perfluor/Fall for Materia Only Pan Perfluor/Fall for Materia Only	40			- 3.0 - 2.5 - 2.0 - 1.5
Intensity Calculations (cm Create Material P 8. A Curreat 8. B Std Correat Time (in seconds per simulations (in o	eate PAR files for one or all a like in PENEPMA_Path	AT files) ("10 hours each at 3600 s 80wwe	40 40			- 3.0 - 2.5 - 2.0 - 1.5
Internity Calculations (cr. Create Material F 8. 8 Commit 8. 8 Statement Time (in seconds per simul Number of Simulations (in 7. Automatically Adjust Electron/Phot	exter PAR files for one or all a like in PENEPHA_Path	AT files) ("10 hours each at 3500 s Bowe at the second state of th	40 40			- 3.0 - 2.5 - 2.0 - 1.5 - 1.0
many Internity Calculations (cre Create Material P R. A Current B Stall Corrent R. B Stall Corrent Time (in seconds per simul Number of Simulations (in of Meinzum Dieston Energy (RM)	elle : PAB files for one or all .8 liles in PENEPHA_Path ellon, x10)	AT Tites) ("10 hours each at 3500 s Bowe AL Boes mathical AT AL Boes mathical Bowe Barren AL Boes mathical Bowe Barren And Al	4 A			- 3.0 - 2.5 - 2.0 - 1.5 - 1.0
many Intensity Calculations (cm Crede Material F Cannot R. B Connot Connot Connot Thre (in second per simulations (in of Automatical) Advant Electron/The Minimum Electron Energy (Ref)	ele .PAB files for cen or all ,8 lies in PDNEPMA_Pah alion, x10) 5000 electrons, x100 [2.0e+09 on Minimum Energy for Z < 11 [Na] [0.4	Art Files) (**10 hours each at 3600 st Bowes Bowes Bowes Bowes Bowes Bowes Par Perhozoffal for Materia Orfer Par Perhozoffal for Materia Std Orig	40 40 40			-3.0 -2.5 -2.0 -1.5 -1.0
many Internity Calculations (cre Create Material P R.A. Cramat B.B. Commat R.B.Stal Commat Time (in seconds per simul Number of Simulations (in r Automaticular Adust Electron/Phost Minimum Electron Energy (RM)	ele PAB Riss for one or all a lies in PENEPHA_Path aton, x10)	AAT files) ("10 hours each at 3600 st Bowre Barner AL Ubes and Faal AL Absorbation (generation PAR Here) Bowre Barner Anner Al Al Absorbation PAR Here) Bowre Barner Al Al Absorbation PAR Here) Barner Perkonfrad for Materia Stat Only Incerner Al Al Absorbation States and Al	δα δΑ δ0 40 40 40 40 40 40 40 40 40 4			- 3.0 - 2.5 - 2.0 - 1.5 - 1.0 - 0.5
many Intensity Calculations (cre Control Matrix) C 4. A Current 4. B SM (Consult Time (n seconds per simul Number of Simulations (in Minisum Electron/Print Minisum Electron/Energy (keV)	techter, x10) [050] techter, x10) [050] techterst, x10) [20x-09] on Minimut Energy to Z < 11 [No] [0.4] K-ratio Calculations: Couple	AT files] ("To hours each at 3000 s Bowers) ("Ben Penkurs end Fran 1 Aff. Ideas materials Bowers) ("Ben Penkur fran 1 Pan Penkur fran 1 Pa	for 3 Å 4 Å 4 Å 4 Å 4 Å 4 Å 4 Å 4 Å 4			- 3.0 - 2.5 - 2.0 - 1.5 - 1.0 - 0.5

Further corrections for Bragg defocussing effects which can vary from negligible to significant depending on the analyzing crystal resolution and orientation of the WDS spectrometer and direction of the fluorescing phase boundary are currently being implemented.

These secondary boundary fluorescence artifacts can also be modeled and corrected "off-line", using our new Penepma 2012 GUI feature in CalcZAF which can be downloaded for free from our web site.

http://www.probesoftware.com/Technical.htm